

## DATASHEET

### Pyrene Actin (10%)

From rabbit skeletal muscle  
( $\alpha$ -skeletal muscle actin, >99% purity)

For Use in Research Only.  
Not for Use in Diagnostic Processes.

Quantity: 1x1.0mg / 5x1.0mg / 10x1.0mg  
Cat. #: 8102-01

#### Product Description

G-actin from rabbit skeletal muscle is a single chain polypeptide with a molecular weight of 42kD consisting of 375 amino acids. Pyrene Actin is chemically modified G-actin carrying a pyrenyl group covalently bound to Cys374 at the C-term of the molecule. Pyrene Actin (10%) contains 10% of pyrenyl actin with a dye/protein molar ratio of 1.0.

The product is supplied as a lyophilized powder containing 2mM Tris-Cl pH 8.2, 0.4mM ATP, 0.5mM DTT, 0.1mM CaCl<sub>2</sub>, 1mM NaN<sub>3</sub> and 0.3% disaccharides, when reconstituted with 1ml H<sub>2</sub>O. For dilution of Pyrene G-actin or exchange of ligand buffer into an actin compatible buffer, MonoMix (Cat.#: 5100-01) may be used. In the absence of nucleators, the polymerization of G-actin can be initiated by PolyMix (Cat.#: 5000-01).

Due to a proprietary cryoconservation technique this highly soluble Pyrene Actin is of excellent biological activity and quality.

#### Product Handling

##### Preparation of a working stock

Example: add 1ml of H<sub>2</sub>O to the tube with 1mg Pyrene Actin to obtain a stock solution of 1mg/ml (24 $\mu$ M). A stock concentration of 4mg/ml (96 $\mu$ M) should not be exceeded to avoid spontaneous polymerization. Allow the G-actin solution to rehydrate for 2min at room temperature and dissolve using a pipette. If Pyrene Actin is used for polymerization with PolyMix (Cat.#: 5000-0\*), we recommend to prepare a G-actin stock of 1.1mg/ml (26 $\mu$ M), as addition of PolyMix will dilute the G-actin/F-actin stock to a final concentration of 1mg/ml (24 $\mu$ M).

For fluorometrical measurements Pyrene Actin may be dialyzed o/n against 50-100vol. of MonoMix or a G-actin compatible buffer and subsequently centrifuged at 100.000xg for 60min to remove any aggregates from freeze-drying.

*For product inquiries please contact:*

cusserv@hypermol.com  
Fon: +49 (0)521 9876228  
Fax: +49 (0)521 9876231  
www.hypermol.com

## DATASHEET

### Pyrene Actin (10%)

From rabbit skeletal muscle  
( $\alpha$ -skeletal muscle actin, >99% purity)

For Use in Research Only.  
Not for Use in Diagnostic Processes.

#### Product Storage and Stability

Store this product at  $-70^{\circ}$  upon arrival, where it will be stable for at least 6 months. Once dissolved Pyrene Actin is kept on ice protected from light and will be stable for at least 4 days. Avoid refreezing of Pyrene Actin solutions.

#### Selection of Related Products

<u>Product</u>	<u>Quantity</u>	<u>Cat. #</u>
Actin (alpha-Actin skeletal muscle, rabbit)	4x250 $\mu$ g	8101-01
Actin (alpha-Actin skeletal muscle, rabbit)	1x1.0mg	8101-03
Atto488-Actin (alpha-Actin skeletal muscle, rabbit)	1x100 $\mu$ g	8153-01
Atto488-Actin (alpha-Actin skeletal muscle, rabbit)	5x100 $\mu$ g	8153-02
Pyrene-Actin 100% (alpha-Actin skeletal muscle, rabbit)	1x1.0mg	8112-01
PolyMix (10x) F-actin buffer	1x1.0ml	5000-01
ME-Buffer (10x) Magnesium Exchange Buffer	2x1.0ml	5111-01
KMEI-Buffer (10x) Magnesium Exchange Buffer	2x1.0ml	5030-01

More products and further information on [www.hypermol.com](http://www.hypermol.com).  
We welcome your visit!

HYPERMOL EK WARRANTS THAT ITS PRODUCT ARE CONFORM TO THE INFORMATION CONTAINED IN THIS PUBLICATION. THE PURCHASERS MUST DETERMINE THE SUITABILITY OF THE PRODUCT(S) FOR THEIR PARTICULAR USE. ADDITIONAL TERMS AND CONDITIONS MAY APPLY (SEE REVERSE SIDE OF THE INVOICE). HYPERMOL BRANDS ARE SOLD EXCLUSIVELY BY HYPERMOL EK, GERMANY AND AUTHORIZED DISTRIBUTORS.

*For product inquiries please contact:*

cusserv@hypermol.com  
Fon: +49 (0)521 9876228  
Fax: +49 (0)521 9876231  
[www.hypermol.com](http://www.hypermol.com)